

## **60° DM Instructions**

Up-dated September 2013

### Personal Protective Equipment:

- a. Lab Coat
- b. Safety glasses/goggles

1. Weigh and record pan # and weight
2. Tare scale
3. Weigh ~100g of sample into pan
4. Record weight
5. Create a duplicate of sample
6. Put in oven for 48 hours
7. Remove and weigh samples back
8. Record weights

Samples to save WET (1 whirl pak/sample)

SFC Sorghum Silage

HMC Corn Silage

WCGF

Samples to save DRY (1 whirl pak/sample both pans)

ALL Mead Feed samples